

hyperClone® Human Fresh Blood
NK Activation/Expansion Kit

Kit05



人新鲜血样来源NK增强版激活扩增试剂盒

NK > 产品性能



新鲜血样来源NK细胞，不受外周血和脐带血起始样本中NK细胞纯度的个体差异性的影响，纯度稳定 > 90%，确保NK细胞培养成功率和质量放行达标

外周血NK以自体使用为主，可以允许有一定量（< 10%）的T细胞存在



大健康NK制剂，追求高性价比

一份50ML新鲜血样培养一次供大健康客户使用

NK > 外周血新鲜血样培养

培养时间	体积 (ml)	计数*10 ⁶	总数*10 ⁶	增值倍数	操作
D0	10	0.39	3.91	1	T25, 10ml+10%自体血浆
D3	10	-	-	-	T75, 补液20ml+10%自体血浆
D5	30	1.83	54.9	14.04	T175, 补液60ml+5%自体血浆
D6	90	2.15	193.5	49.49	T175,补液50ml+5%自体血浆
D7	150	2.23	334.5	85.54	入袋, 补液150ml+自体血浆2%
D8	300	2.46	738	188.75	补液300ml
D9	600	2.12	1272	325.31	补液600ml
D11	1200	2.33	2796	715.09	补液600ml
D14	1800	2.84	5112	1307.41	收获

NK > 数据展示

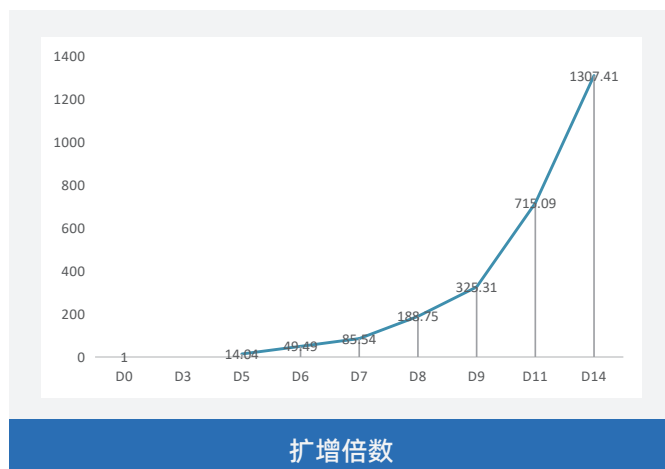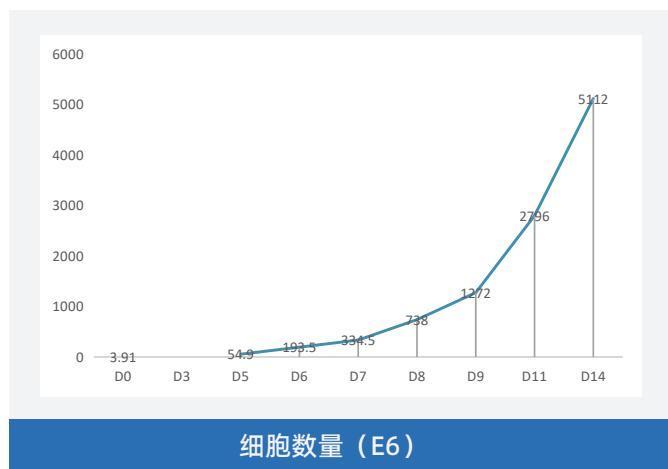

NK > D-14天流式图

